


DRD's patented **Differential Resolution Displacement** pistons give extremely fine resolution over a very wide range along with fast flow and large volume. The 3rd generation, trademarked **CV1dur™**, drives the pistons directly and coaxially with the integral lead screw of the latest and best linear actuator motor technology.

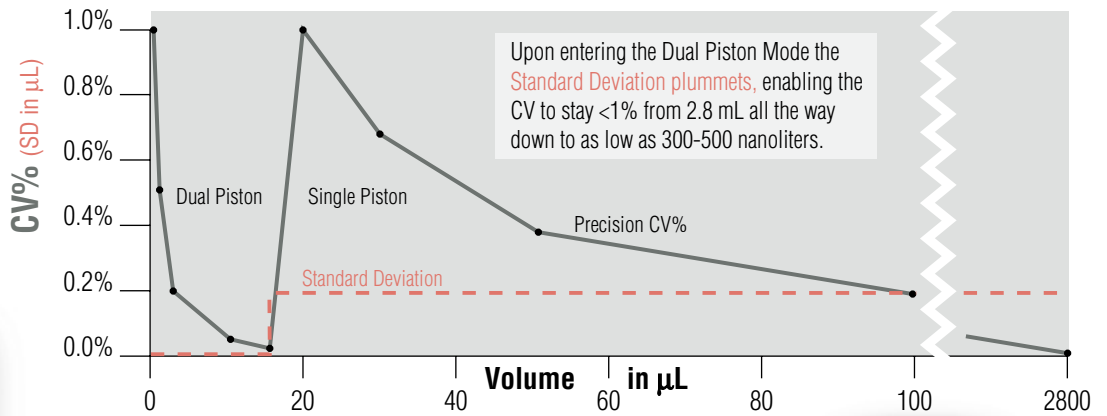


DUAL PISTON MODE

When the spring-loaded top piston, which is of slightly smaller diameter than the bottom piston, follows the bottom piston and **the two pistons move together** in the upper part of the chamber, the volume of liquid metered equals the small difference in the volume displaced by the two pistons (thin annular ring). Volume resolution is extremely fine and gives **high precision**, reflected in CVs that hold below 1% to extremely low volumes.



SINGLE PISTON MODE

When the motor-controlled **bottom piston moves alone** in the bottom part of the chamber, the volume of liquid metered equals the volume displaced by only the bottom piston (solid wafer), giving **large volume** and **fast flow** rate. This provides easy priming, bubble clearance, clean shear-off, and discrete tiny sample pick-up and touchless **Blastoff™** delivery.

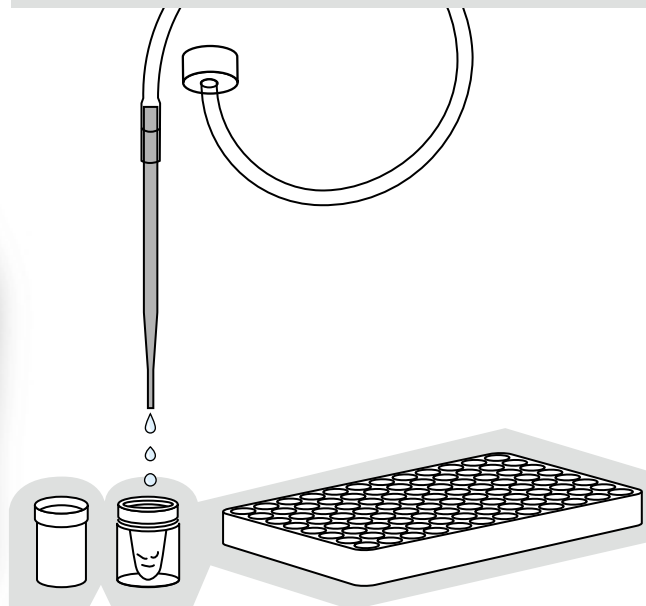

MODEL 48T ACTUAL SIZE
8.9" (22.6 cm) high



MODEL 48S ACTUAL SIZE
6.1" (15.5 cm) high