


# DRD NanoBlast96™

Pipetting full range with Model Anthony and Labcon50 disposable tip

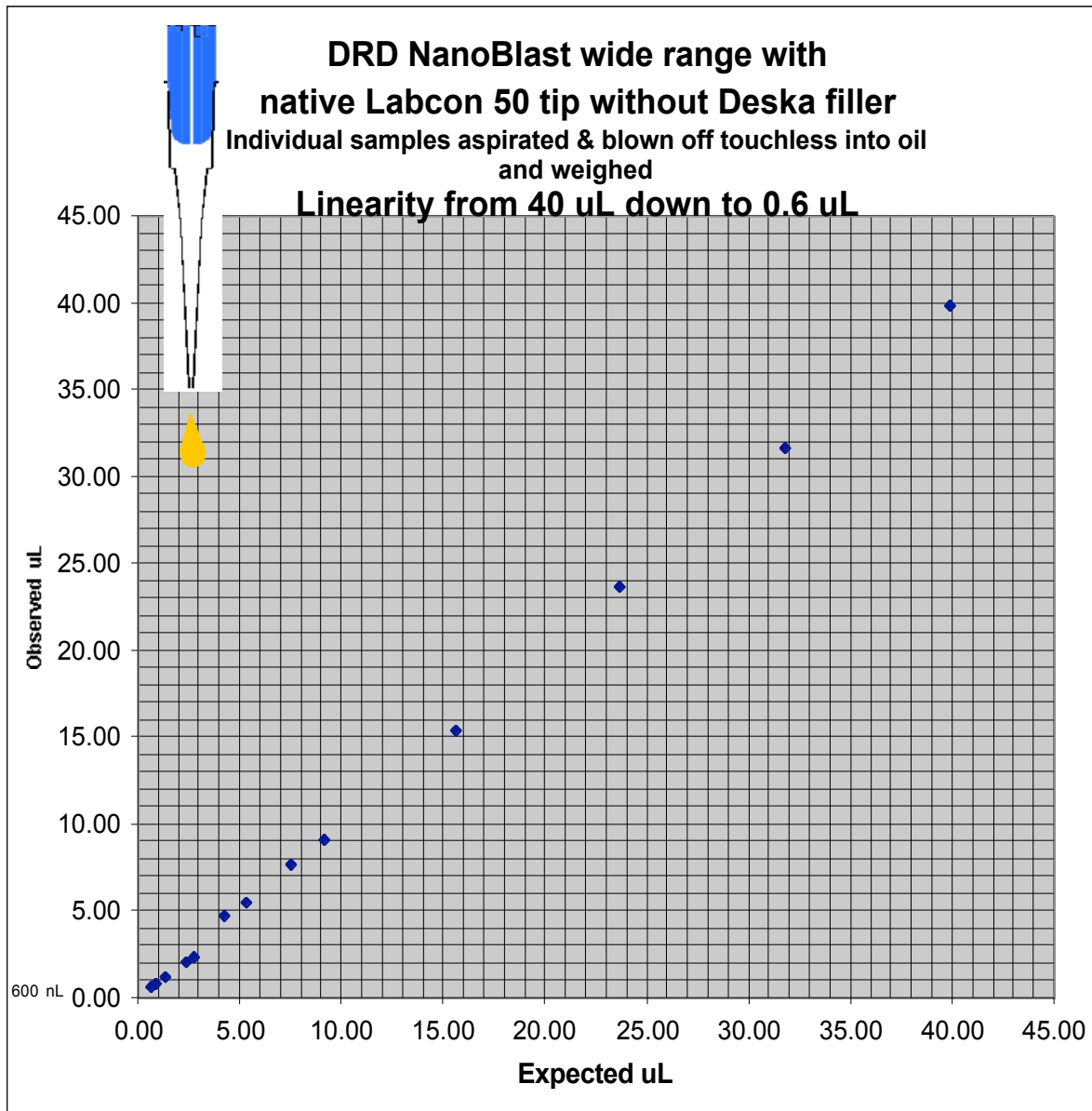

The samples over 5 uL were aspirated in the Bulk or Single Piston mode and those below 5 uL were aspirated in the fine Differential Mode. After aspiration of each discrete sample, the sample was blown off in the Bulk mode -- contact free -- into a cup of oil and weighed.

The linear relationship extends down to 500 - 600 nanoliters.

To see the discrete pipetting to lower nanoliter volumes, see companion "Pipetting nano range and lower limits using Deska filler".